

Fetal Bovine Serum

CAT N°: FB-1001, FB-1003, FB-1058, FB-1061, FB-1280, FB-1285, FB-1290, FB-1295, FB-1345, FB-1350, FB-1360, FB-1365

Collected from the source:

When searchers choose their serum an important factor that should be taken into consideration is the source, which also emphasises the traceability of the serum.

Our system of vertical integration allows us to be certain of the origins and traceability of our FBS.

Each manufactured batch is rigorously controlled, from the collection of serum and throughout all stages of its treatment and production through to final packaging on our premises.

BioSera Fetal Bovine Serum is derived from clotted whole blood aseptically collected from foetus via cardiac puncture.

The serum is collected or imported and treated in agreement with the European regulations.

Country of Origin:

The country in which the serum was taken from the donor/animal.

<u>Cat. N°</u>	<u>Description</u>	<u>Origin(s)</u>
FB-1001	Fetal Bovine Serum (South America Origin)	South America
FB-1003	Fetal Bovine Serum (South Africa Origin)	South Africa
FB-1058	Fetal Bovine Serum (Uruguay Origin)	Uruguay
FB-1061	Fetal Bovine Serum (Dominican Republic Origin)	Dominican Republic
FB-1280	Fetal Bovine Serum (French Origin)	France
FB-1285	Fetal Bovine Serum (Ireland Origin)	Ireland
FB-1290	Fetal Bovine Serum (Spain Origin)	Spain
FB-1295	Fetal Bovine Serum (Netherland Origin)	Netherlands
FB-1345	Fetal Bovine Serum (Central America, USDA approved)	Costa Rica, Panama, Honduras, Guatemala
FB-1350	Fetal Bovine Serum (USA Origin)	United States of America
FB-1360	Fetal Bovine Serum (Mexico, USDA approved)	Mexico
FB-1365	Fetal Bovine Serum (Chile, USDA approved)	Chile

Storage conditions:

- 18°C to - 40°C, protected from light.

Bottles can be stored between -40°C and -80°C for a short period (few days).

Shelf life:

5 years

Filtration:

Final Filter Size: 0.1µm x 3

pH:

pH specification: 7,4 ± 0,6

Osmolality:

Determined by a lowered freezing temperature. The osmometer is calibrated against standard solutions.

Osmolality specification: 322,5 ± 42,5 mOsm/kg

Endotoxin:

All sera are tested to determine the levels of endotoxins. BioSera carries out a chromokinetic quantitative test, according to the method D of the European Pharmacopoeia.

The endotoxin reagent is standardized against the US reference endotoxin.

Specifications available on Certificate of Analysis

Haemoglobin:

The haemoglobin level is measured by spectrophotometer.

Specifications available on Certificate of Analysis

Cell Culture:

Biological performance is assessed using cell culture medium supplemented with the serum being tested.

During the test period, cultures are examined microscopically for any morphological abnormalities that may indicate toxic components in the serum.

Cell Culture Tests:

Cell Growth, Plating Efficiency, Cloning Efficiency.

Cell Lines Tested:

The following cell lines are tested with the serum:

HELA -Cancer Cell/Human.

L929 -Fibroblast-Mouse/ As Macrophage

SP2/0-AG14 -Mouse/Lymphoma

MRC- 5 -Human/Lung.

Total Protein:

Determined by Biuret Colorimetry.

Total Protein specification: 40 ± 15 g/l

Sterility tests:

All sera are tested for the absence of aerobic and anaerobic bacteria, fungi, yeast and *Mycoplasma*.

The sterility test is based on the European Pharmacopoeia requirements.

The sera are tested for the absence of *Mycoplasma* by culture.

Virus test:

All of our sera are tested for:

- Bovine Viral Diarrhoea (BVD)
- Cytopathogenic agents e.g. Infectious Bovine Rhinotracheitis (IBR) / BHV-1
- Hemadsorbing agents e.g. Parainfluenza Type 3 (PI3)

Sera are tested by inoculation to permissive cells. The revelation is made by immunofluorescence for pestiviruses. Cytopathogenic agents and hemadsorbing agents are detected by microscopic observations.

Other tests:

Not Applicable

Treatments:

Not Applicable

Recommended use:

- Respect storage conditions of the serum
- Do not use the serum after its expiry date
- Store serum in an area protected from light
- Manipulate serum in aseptic conditions (e.g.: under laminar air flow)
- Wear clothes adapted to the manipulation of serum to avoid contamination (e.g.: gloves, mask, hygiene cap, overall...)
- In order to preserve all serum qualities, it is recommended to thaw out the flask, to aliquote, then to re-freeze the produced flasks rather than to thaw out and re-freeze the flask at each use.
- It is recommended to use the serum immediately after its thaw out. However, if it is not useful, it is possible to store thaw out serum, at +2°C / +8°C, until 26 weeks without significant decrease of its performances in cell culture.

The product is intended to be used in vitro for research or further manufacturing only and not for use as an Active Pharmaceutical Ingredient or food or animal feed.

Remarks:

The raw serum may be treated (Heat Inactivated, Gamma Irradiated, pH modified) before filtration for different reasons:

- Importation regulation
- Exportation necessity
- Technical or quality aspects

To be informed if your batch is concerned by treatment before filtration,
please contact BioSera.